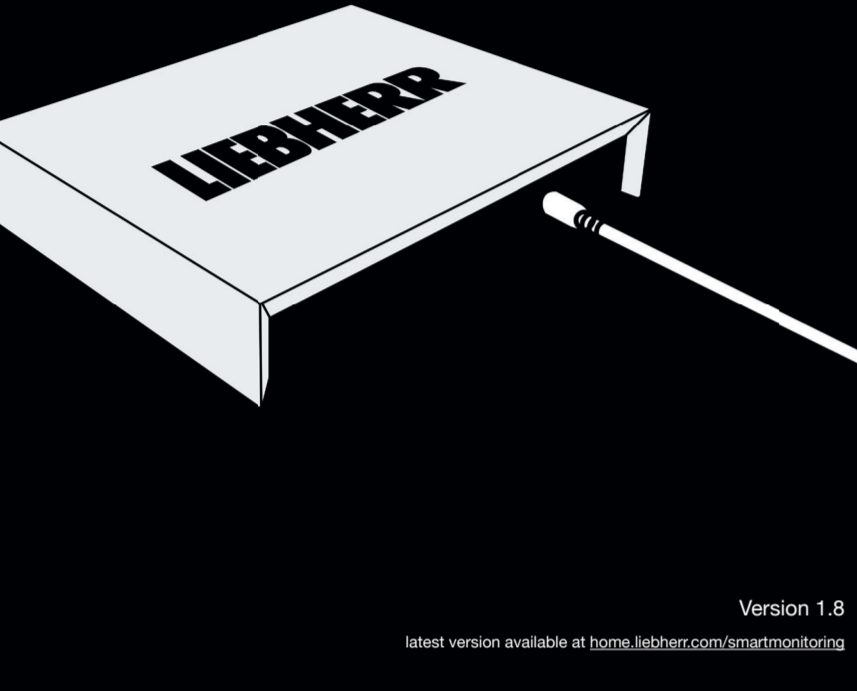
# SmartCoolingHub

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Version 1.8

latest version available at [home.liebherr.com/smartmonitoring](http://home.liebherr.com/smartmonitoring)

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For more information and other accessories visit the SmartMonitoring website: [home.liebherr.com/smartmonitoring](http://home.liebherr.com/smartmonitoring) or contact your dealer.

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<b>Start the set-up at the following address:</b> <a href="http://smartmonitoring.liebherr.com/setup">smartmonitoring.liebherr.com/setup</a> or scan this QR-Code:

You will also find the following there:

- Set-up videos
- Appliance instruction manual
- SmartMonitoring instruction manual

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For more information, refer to the SmartMonitoring instruction manual.

- DHCP mode active (automatic IP address assignment)
- The MAC address of the SmartCoolingHub must be shared in the network (see label on the back of the SmartCoolingHub if a MAC filter is used in your network).
- An Internet-enabled device with a current browser version and a minimum resolution of 1024x768 pixels is required to access SmartMonitoring.
- For your convenience, make a note of the serial numbers of your refrigerators in advance.
- A MyLiebherr customer profile is required to log in to the SmartMonitoring dashboard (registration possible during set-up).
- You will need additional accessories to set-up other appliances.
- Recommendation: one SmartCoolingHub per room
- If the distance between the SmartCoolingHub and the appliance is greater than 50 m, you should consult a qualified technician for installation.

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